


HOW TO USE THE BODE SecurSwab™ FOR BIOLOGICAL STAINS

CATALOG NUMBER P08D71


FOR HUMAN/FORENSIC IDENTIFICATION ONLY—NOT FOR CLINICAL OR DIAGNOSTIC USE

NOTE: Use gloves to prevent sample contamination.

Do not use Bode SecurSwab™ if the Indicating Desiccant pouch is pink.

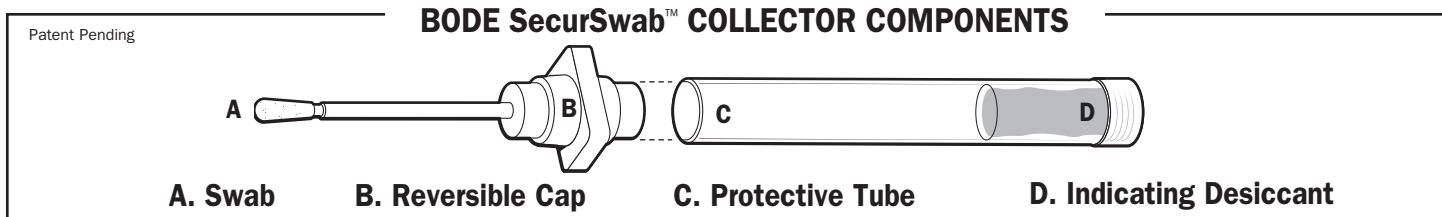


Step 1

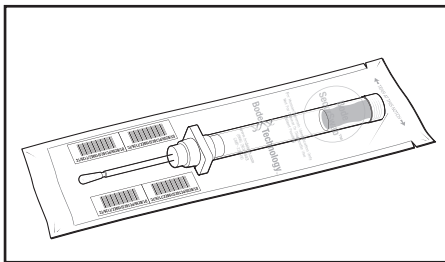


- Place the wrapped collector near the evidence and ensure the barcode label is visible, if photo documentation is required. Make sure that the unique barcode label on the collector package is visible through the camera eye (optional).
- Tear open the collector wrapper at the notches and remove the collector by holding the Protective Tube. Remove one barcode label from the wrapper and apply to the Protective Tube for sample ID. Save the other barcode labels for labeling a transport pouch or appropriate documentation.

Step 2

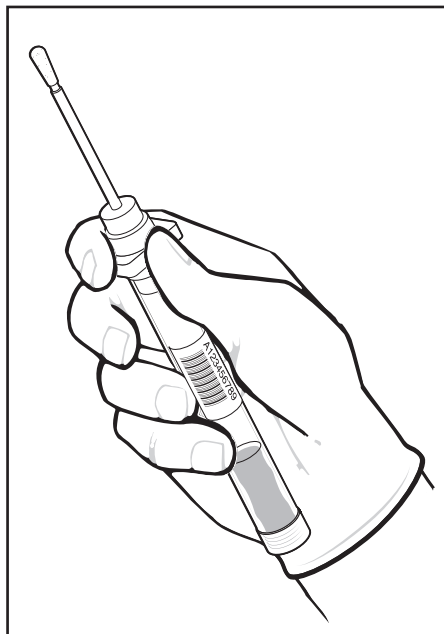


- Hold the collector firmly in your hand, as shown.
- Keep your forefinger over the Reversible Cap, as you collect the sample.

Step 3

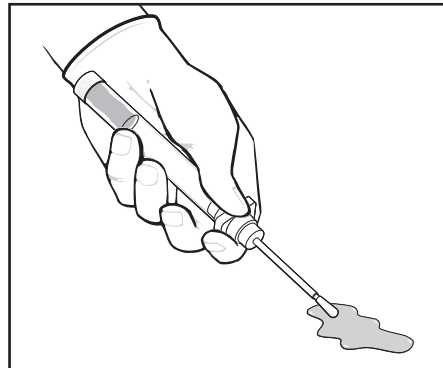


- For dry stain collection: Add 100 microliters of sterile water (about 3 drops) to the swab and collect by touching and rotating the swab on stain.
- For wet evidence collection: Dab the swab in the evidence to collect about 100 microliters (about 3 drops) of evidence. Air drying will be necessary if too much fluid is collected.

Step 4

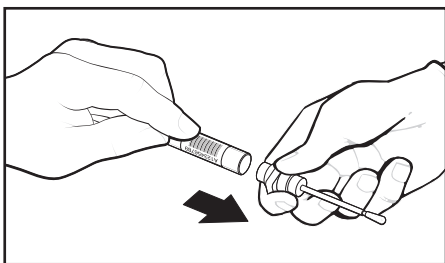


- Firmly grip the Reversible Cap with one hand and the Protective Tube in the other hand.
- Twist and pull the Protective Tube away from the Reversible Cap: Be careful not to touch the Swab.

Step 5

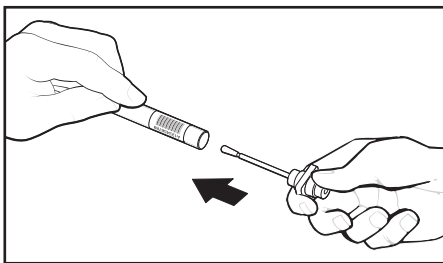

- Carefully turn the Reversible Cap around 180 degrees.
- Insert the Swab in the Protective Tube without touching the Swab to the inside of the Protective Tube.
- Press and twist the Protective Tube firmly into the Reversible Cap to seal.

Step 6

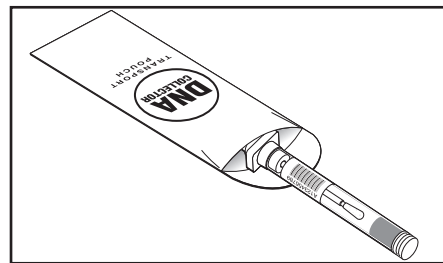

- Insert the collector into a transport pouch and seal.
- Remove barcode labels from the collector wrapper and apply to transport pouch for evidence identification.
- Discard the collector wrapper.


HOW TO USE THE BODE SecurSwab™ FOR BUCCAL CELL COLLECTION

CATALOG NUMBER P08D71


FOR HUMAN/FORENSIC IDENTIFICATION ONLY—NOT FOR CLINICAL OR DIAGNOSTIC USE

NOTE: Use gloves to prevent sample contamination. If you are self-collecting, you are the subject in steps 2–5.

Do not use Bode SecurSwab™, if the Indicating Desiccant pouch is pink.

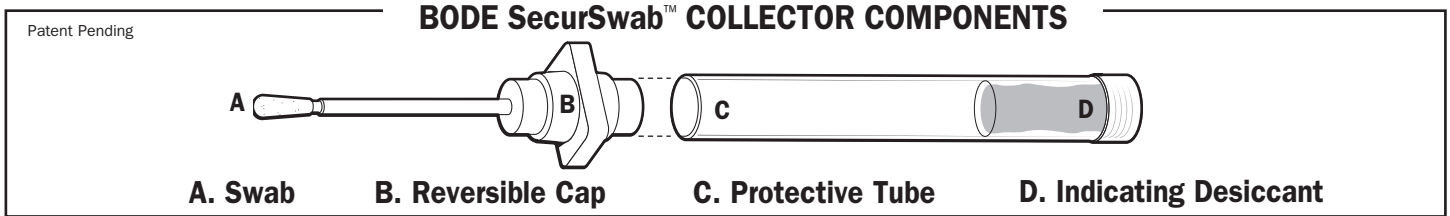


Step 1

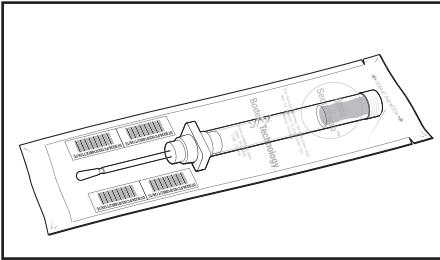

- Tear open the collector wrapper at the notches and remove the collector holding by the Protective Tube.
- Remove one barcode label from the wrapper and apply to the Protective Tube. Save the other barcode labels for labeling a transport pouch or appropriate documentation.

Step 2

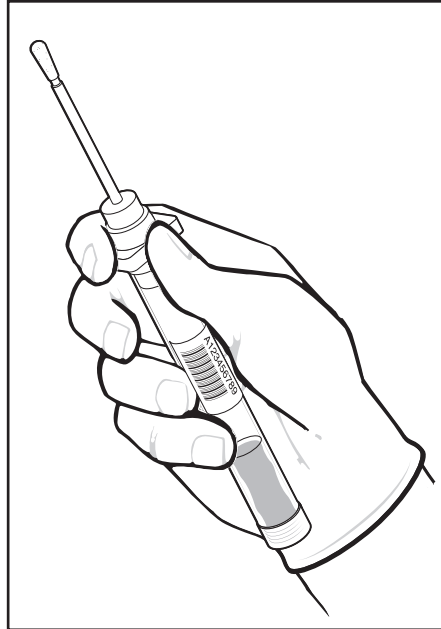


- Hold the collector firmly in your hand, as shown.
- Keep your forefinger over the Reversible Cap, as you collect the sample.

Step 3




- Instruct subject to open mouth and place the swab inside the mouth against the cheek.
- Ask the subject to press the Swab against the cheek and move the Swab in 10 circular patterns, about 1 inch in diameter. At the same time the subject must rotate the Protective Tube to moisten the complete Swab.
- It is important for the collection personnel to observe a bulge in the cheek, as the subject presses and rotates the Swab against the cheek.

Step 4

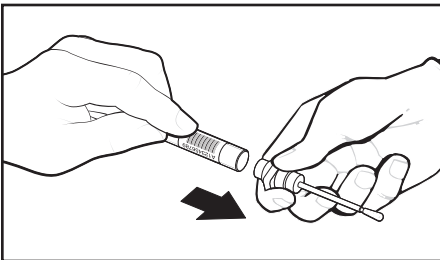


- Firmly grip the Reversible Cap with one hand and the Protective Tube in the other hand.
- Twist and pull the Protective Tube away from the Reversible Cap: Be careful not to touch the Swab.

Step 5

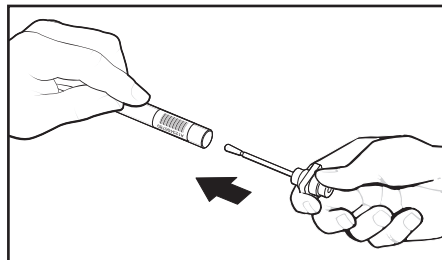

- Carefully turn the Reversible Cap around 180 degrees.
- Insert the Swab in the Protective Tube without touching the Swab to the inside of the Protective Tube.
- Press and twist the Protective Tube firmly into the Reversible Cap to seal.

Step 6

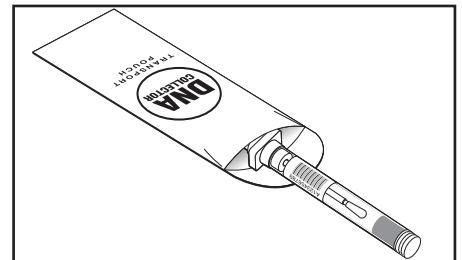

- Insert the collector into a transport pouch and seal.
- Remove barcode labels from the collector wrapper and apply to transport pouch for identification.
- Discard the collector wrapper.