

Bode Buccal 2 Collector

Instructions for Sample Conversion into Cassette Format



Bode Buccal 2 Cassette


A. Clean Punch Area

B. Pegs

C. Guide Rails

D. Bode Buccal 2 Coin Envelope

NOTE: Please allow the buccal sample to completely dry prior to conversion.

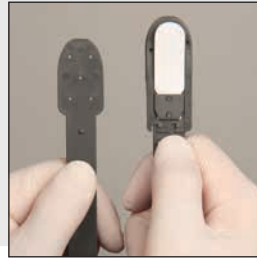

Step 1

The individual converting the collector into the cassette format should wear protective gloves. Fully expose the collection paper by moving the slider downward. Remove the slider from the Bode Buccal 2 by pulling it apart from the handle. The slider can be discarded according to the laboratory protocol.



Step 2

Remove the Bode Buccal 2 cassette from the coin envelope (D). Place the cassette in one hand while holding the Bode Buccal 2 in the opposite hand with the THUMB designation facing you.



Step 3

Align the bottom left and the bottom right edge of collector frame with guide rails (C) on the cassette. Slide the collector frame until it stops.



Step 4

Place the cassette/collector on a clean surface. Press the top right edge of the handle down until the peg clicks into the collector frame. Repeat this action on the top left edge.



Step 5

Slightly lift the front end of the handle, creating a gap between the handle and the collector frame. Applying pressure on the grips, shift the handle to the right and pull down to disconnect the handle.

Note: If the pegs are not pushed completely through the collector frame holes, the edge of the handle can be used for assistance.



Step 6

Discard the handle according to the laboratory protocol and proceed with sample processing or storage in the desiccated foil pouch.